


**Perfusion Pressure Kit (#09-04, 09-08 & 09-16)**

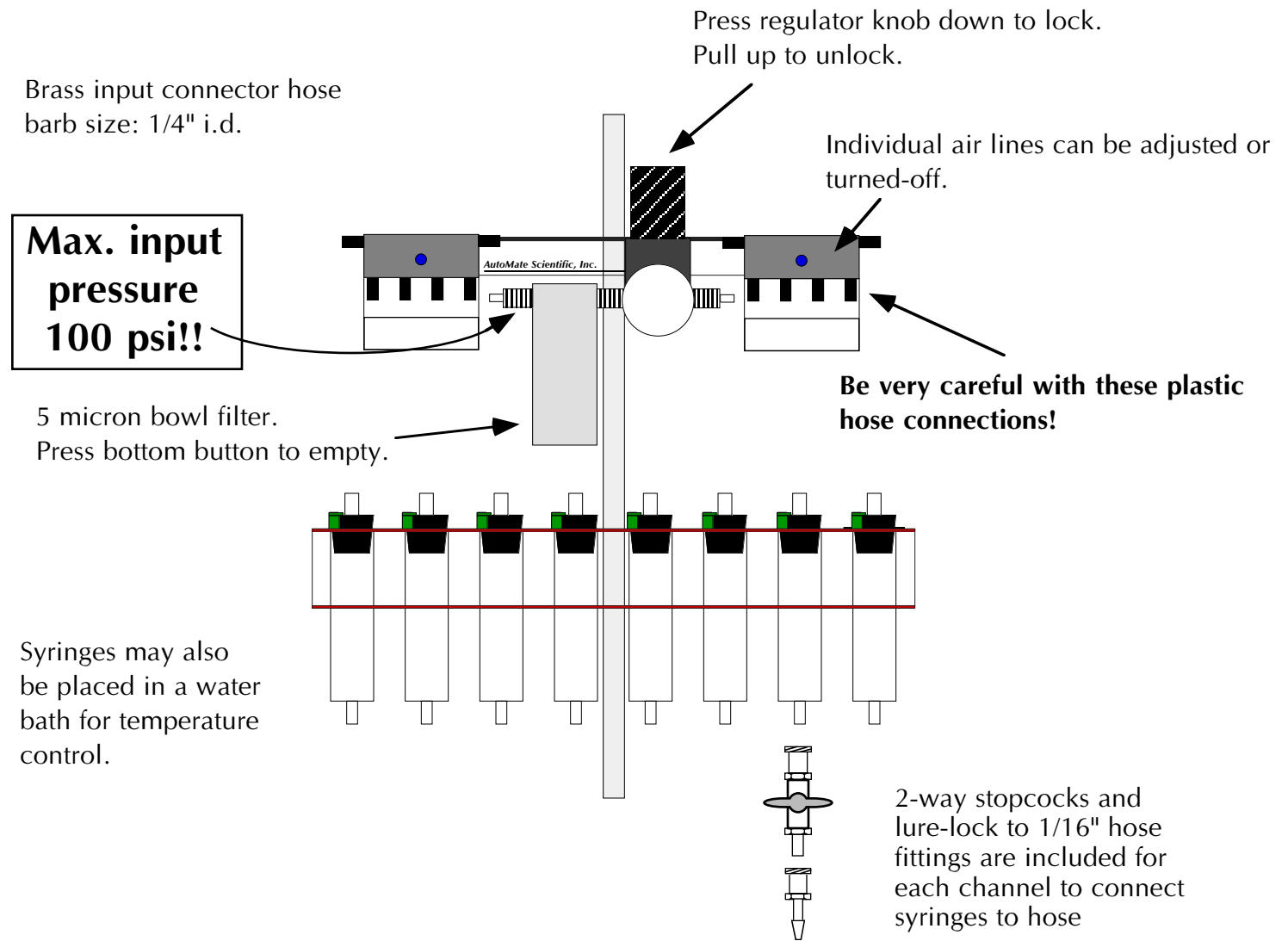

Brass input connector hose  
barb size: 1/4" i.d.

Press regulator knob down to lock.  
Pull up to unlock.

**Max. input  
pressure  
100 psi!!**

Individual air lines can be adjusted or  
turned-off.

**Be very careful with these plastic  
hose connections!**

5 micron bowl filter.  
Press bottom button to empty.

Syringes may also  
be placed in a water  
bath for temperature  
control.

2-way stopcocks and  
lure-lock to 1/16" hose  
fittings are included for  
each channel to connect  
syringes to hose

Hoses should continue from syringes to valves.



800.998.MATE | [www.autom8.com](http://www.autom8.com) | 650 University Ave #5, Berkeley, CA 94710 USA  
tel 510.845.6283 | fax 510.665.3975 | e-mail [info@autom8.com](mailto:info@autom8.com)